


Acquire and Analyze Q-Plex Arrays

The Q-View Software is a powerful tool for the high sensitivity imaging and quantitative analysis of Q-Plex Arrays and other assays, such as classic ELISAs and blots. It drives the Q-View Imager, applying advanced image acquisition algorithms to minimize the effects of noise, maximize sensitivity, and eliminate the need for exposure stacking. Q-View Software offers many time-saving "Auto" features, multiple regression models including 4PL and 5PL, three weighting options to offset heteroscedasticity, outlier masking, pass/fail calculations to automatically detect and flag errors, and customizable reports and charts. Related images and their large data sets are housed within a single project file and exported for further use.

Intuitive Workflow

The software enables data analysis in three easy steps:

- 1. Image Processing:** Acquire or import images and easily locate spots using Auto-Set Plate Overlay and Auto-Adjust Spots Or use Blot Tools to determine signal intensities from alternate assay types.
- 2. Well Assignment:** Label wells quickly via Templates and Sequential Naming, and import/export plate layouts as .csv files.
- 3. Data Analysis:** View raw data, Auto-Select the best curve fit or manually optimize regression model settings for your data, and customize charts and reports with Statistics, Assay Limits, and Auto-Selected Dilutions.

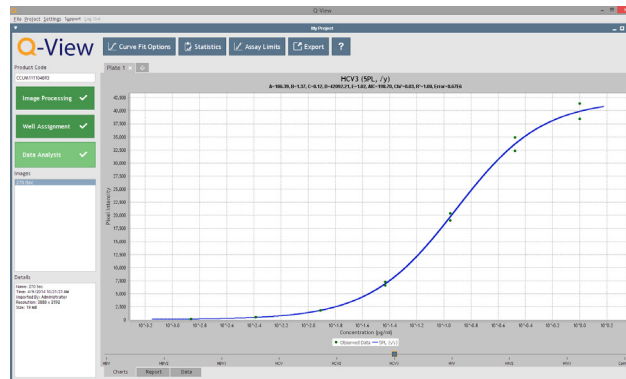

Advantages

- ▶ cGMP-compatible
- ▶ Full touch-screen support
- ▶ Concise embedded help and online video tutorials
- ▶ Publication-quality, customizable reports and graphs
- ▶ Rapid processing through multi-threading and task parallelization

Computer Requirements

	Minimum	Recommended
Operating System: (32 or 64 bit)	Windows 10, 8.1, 8, 7, Vista or XP	Windows 10, 8.1, 8, 7, Vista or XP
Ram:	2 GB	3+ GB
Java:	32-bit Version 7 or greater	Current 32-bit Version
Processor:		2 or more processing cores

Compatible Instruments

The following chemiluminescent instruments are engineered for maximum optical performance with Q-Plex Arrays and automated via Q-View Software.

- ▶ Q-View Imager Pro
- ▶ Q-View Imager LS

How to Get the Q-View Software

Q-View Software can be installed on any number of computers. To request a quote for Q-View Software, contact your Quansys Sales representative (888-782-6797), email info@quansysbio.com, or go online to www.quansysbio.com/q-view-software.



Quansys Biosciences
365 North 500 West
Logan, UT 84321
www.quansysbio.com

Fax: 435-750-6869
Phone: 435-752-0531
Toll Free: 888-QUANSYS (782-6797)

Quansys is an ISO 9001:2008 and ISO 13485:2003 registered company and complies with cGMP. Products are designed, developed, and manufactured according to the procedures outlined in our Quality Management System.

©2018 Quansys Biosciences Inc. All rights reserved. Q-View and Q-Plex are Trademarks of Quansys Biosciences Inc. ISO is a trademark of the International Standards Organization. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

Increase productivity with
Quansys products and services.
Save Time, Sample and Money.

Q-Plex™ ARRAYS

Our Q-Plex Arrays are quantitative multiplex ELISAs with distinct proteins deposited in a defined array. Choose one of our standard kits for immediate delivery or customize the exact array you need.

Q-View™ IMAGERS

The Q-View Imager Pro and Q-View Imager LS are high quality, low-cost imaging systems for chemiluminescent assay imaging.

Q-View™ SOFTWARE

The Q-View Software is a user-friendly image analysis app that enables the acquisition and analysis of large amounts of multiplex ELISA data.

SAMPLE TESTING SERVICE

Our lab runs immunoassays on a wide variety of biological samples. Increase productivity and focus on your core research while we test your samples for you.

QUANSYS
B I O S C I E N C E S



www.quansysbio.com